



Houston Forensic Science Center

Comparative and Analytical Division - Toxicology

Hamilton® Maintenance Log

Manufacturer: Hamilton Company

Model: Microlab® 600

Serial Number: 1742 or 7933

| Type of Maintenance* | Recommended maintenance schedule* | Check if performed |
|--|--|------------------------------|
| Replacing Valves | Change every 1.5 to 2 years. | ✓ |
| Replacing Cross Tube | Change when valves are changed. | ✓ |
| Replacing the Syringes | Change if solution is leaking past the plunger. | Reinstalled gv 08/29/2023 |
| Replace Tubing | When becomes bent/kinked or discolored or with annual service. | ✓ |
| Cleaning and Lube Service | Annually by Hamilton Company Service Center. | |
| Replacing Batteries | Battery should have a lifespan of 10 years. When battery fails send controller to Hamilton Company Service Center. | |
| Updating Firmware or Software Version | As updated by Hamilton Company | |
| | Version | |
| | Firmware FRGA | gv 08/29/2023 |
| | Firmware Runtime | |
| | Software | |
| Operating System | | |
| gv 08/29/2023 | | |
| Comments: Received back from Hamilton Company on 08/28/2023 - gv 08/29/2023 | | |

*Adapted from Hamilton® user manual.

Signature: _____
 Date: 08/29/2023