



Houston Forensic Science Center

Comparative and Analytical Division - Toxicology

Millipore Maintenance Log

Manufacturer: Millipore
Model: Direct-Q 3 UV with Pump
Serial Number: F9NA15328D

Type*	When maintenance is performed*	Initials/date performed
Sanitization of the System	Once a year (SmartPak will be replaced after the sanitization of the system).	/
SmartPak Replacement	1. When the Pack Alarm display is blinking, 2. When the system resistivity display is blinking, or 3. After a system or tank sanitization.	
Tank Level Calibration	When the SmartPak is replaced.	
Screen Filter Cleaning	Recommended 1. When the SmartPak is replaced, or 2. When the filter is clogged.	
Vent Filter Replacement	1. When the SmartPak is replaced, or 2. After a system or tank sanitization.	BAM 4/7/21
Final Filter Replacement	1. When SmartPak is replaced, 2. When the Product Water flowrate drops, or 3. After a system or tank sanitization.	/
Flow Calibration	When the Final Filter is replaced.	
UV Lamp Replacement and UV Timer Reset	When the UV Lamp Alarm display is blinking.	/
Sanitization of the Tank	Recommended once a year (not needed if sanitization of the system has been performed).	
Comments: Sanitized the tank due to the smart pak replacement alarm still showing. -BAM 4/7/21		

*Refer to user's manual for additional details.

Signature: Brooke Mandenhall

Date: 4/7/21